


User guide


NanoDrop Lite Plus quick start guide

NanoDrop Lite Plus Home Screen

From the Home Screen, select the **Nucleic Acids** or **Protein** sample category, then select a sample type from the available applications

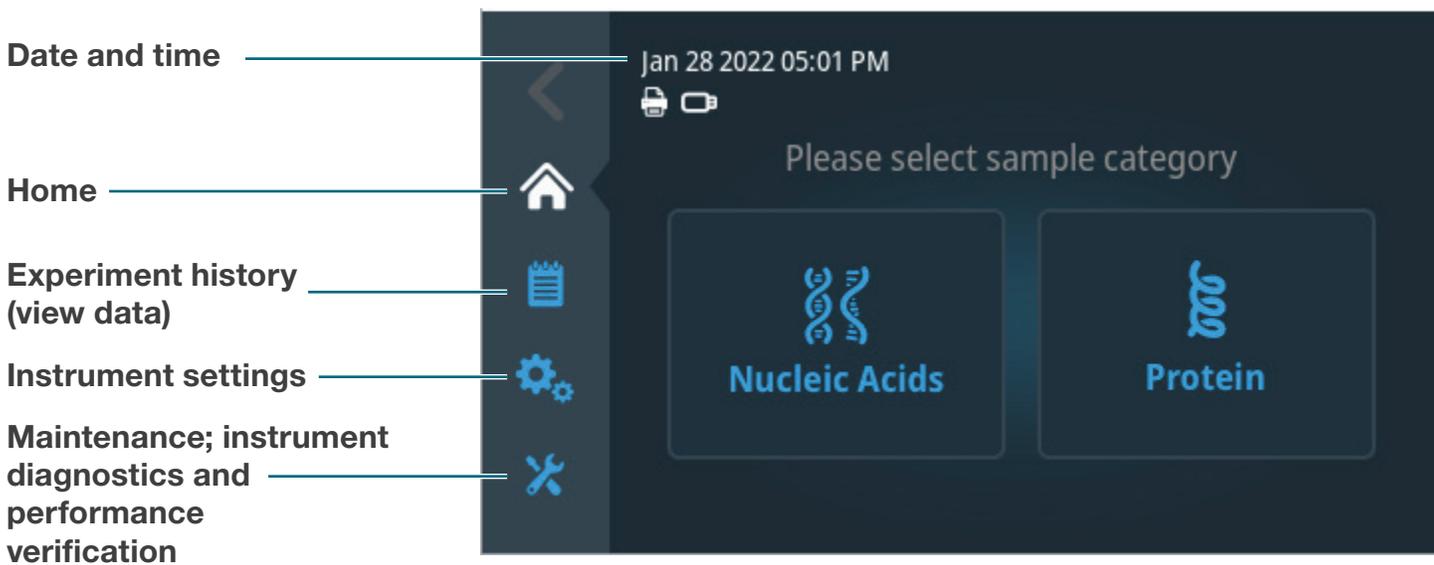


Load blank

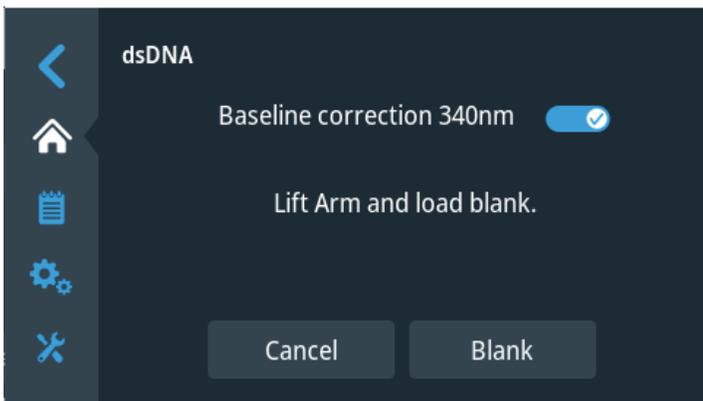


Load blank, close the arm quickly, and select **Blank**.

Run samples

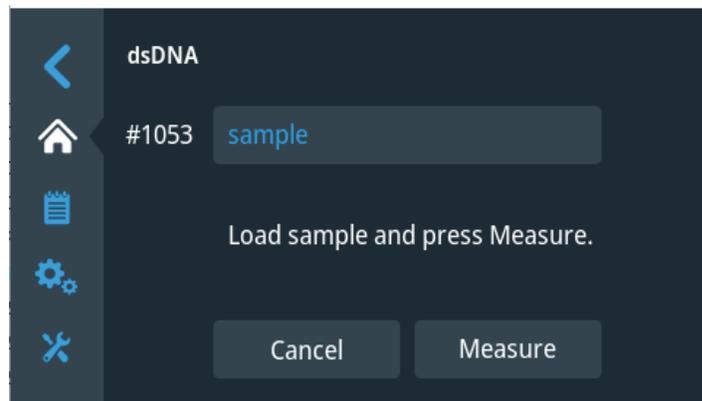

Enter a sample name, load sample, close the arm quickly, and select **Measure** to initiate measurement.

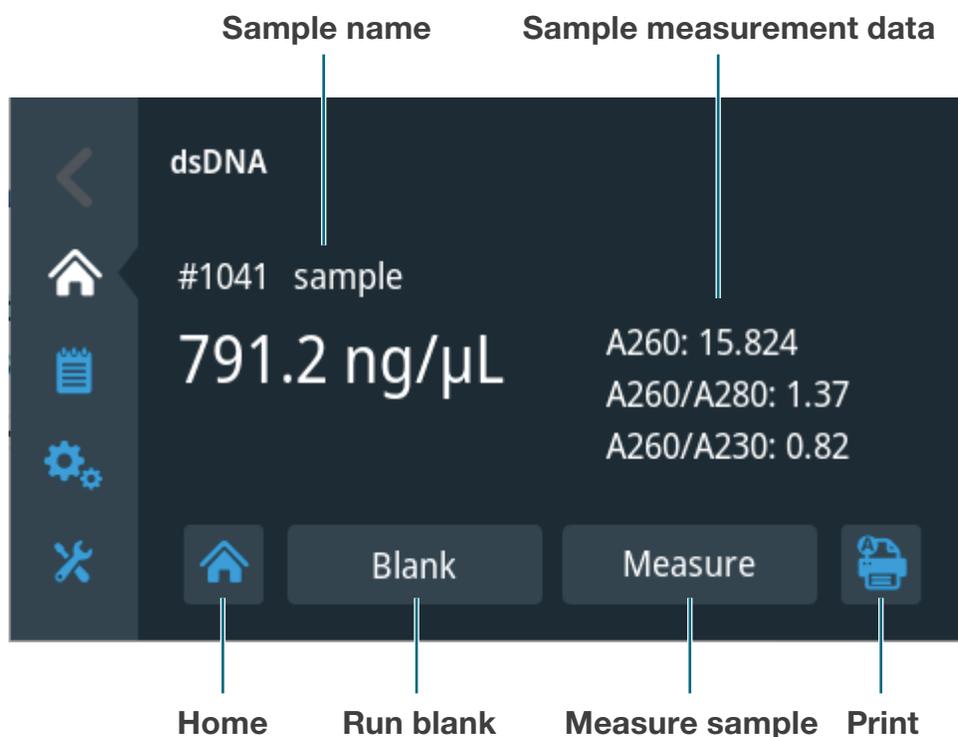

When finished measuring samples, select **Home**. Sample data is automatically saved.

Thermo Scientific NanoDrop Products
302-479-7707

Wilmington, Delaware USA
Toll Free (US & Canada)
877-724-7690

Technical Support:
nanodrop@thermofisher.com
www.thermoscientific.com/nanodrop

Once printed this document is no longer controlled

Learn more at [thermofisher.com/nanodrop](https://www.thermofisher.com/nanodrop)

thermo scientific